


INSTRUCTION SHEET

HOW TO USE GATTAquant Nanorulers


STEP 01

GATTAquant nanorulers are shipped immobilized on a coverslip and are embedded in polymer or imaging buffer.

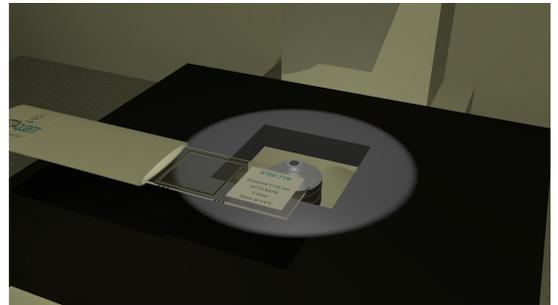

STEP 02

You just put the GATTAquant nanoruler sample on your microscope but make sure that the coverslip side faces the objective (right inset of the picture).

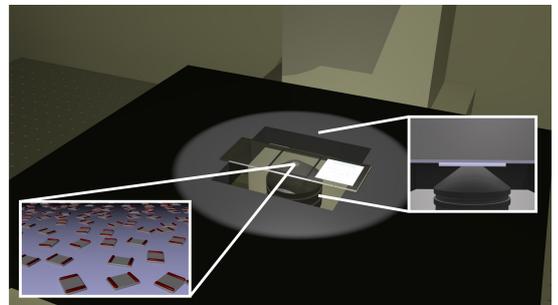

STEP 03

Switch on **laser** and **camera** and **drive carefully the objective control in one direction** until the surface is visible.

We strongly recommend approaching using the **camera** with a rather small field of view ($40\ \mu\text{m} \times 40\ \mu\text{m}$ or smaller). Be aware of that nanorulers are **nanoscaled** and smaller/dimmer than, for instance, stained cells, consequently approaching via the ocular is not advised.

